

IncuSafe FAQs

Hydrogen Peroxide (H₂O₂) Vapor Decontamination



MCO-230AICUVH-PE



MCO-170MUVH-PE




MCO-170AICUVH-PE

H₂O₂ Decontamination Cycle

Hydrogen Peroxide (H₂O₂) Vapor Decontamination is available as an option on these IncuSafe incubators. This system eliminates 100% of contamination. Developed with safe and effective cell culture decontamination processes using H₂O₂ vapor in conjunction with narrow-bandwidth ultraviolet light and copper enriched stainless steel to broaden CO₂ incubator application to highly regulated, environmentally sensitive cell cultures.

Throughout the entire cycle the incubator airflow system continues to gently circulate interior air assuring 100% vapor contact with all interior surfaces, ultimately creating a breakdown of H₂O₂ to harmless water and oxygen as it passes over the UV lamp. Ultimately converting all H₂O₂ to less than 1 part per million following completion of the decontamination cycle. No ozone is produced for this decontamination process.

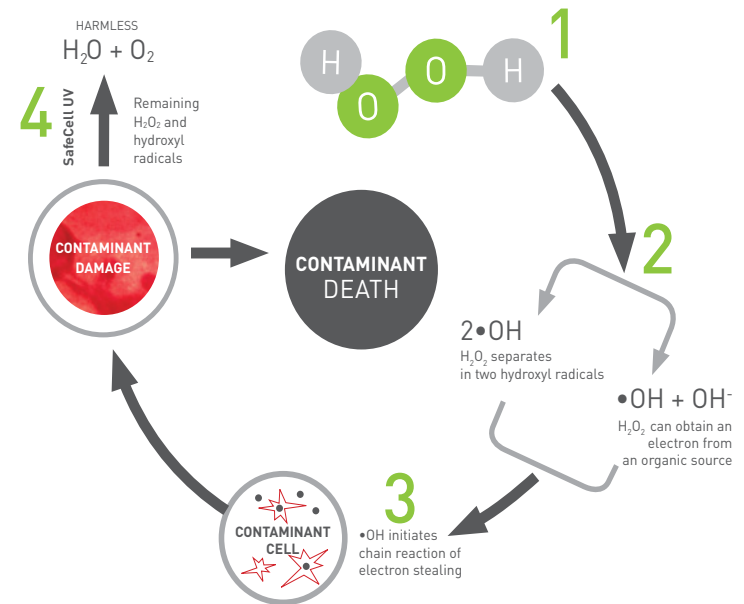

H₂O₂ Section

[What do I need to do if power failure occurs during the decontamination process?](#)

[How is H₂O₂ resolved?](#)

[What marks a resolving phase completion?](#)

[During decontamination and resolving processes, is the chamber fan operating?](#)

[What triggers the switch from H₂O₂ generation to the resolving phase?](#)

[After decontamination, is it OK to leave the incubator unattended for a night or a couple of days? \(Can I start the decontamination program just before going home?\)](#)

[Can I use a commercial H₂O₂ solution?](#)

[When should I perform the decontamination process?](#)

[Why do you use a 6% MCO-H2O2-PE solution?](#)

[What temperature is needed for the storage of the H₂O₂ reagent?](#)

[How should I store H₂O₂ reagent \(MCO-H2O2-PE\)?](#)

[What happens if MCO-H2O2-PE is frozen? Is it OK to put it in a freezer when the expiration date has been exceeded?](#)

[How is H₂O₂ reagent transported?](#)

[Is it possible to put something that I want to decontaminate into the MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE incubators and decontaminate it with H₂O₂?](#)

[How is decontamination affected if extra trays or half trays are added into the incubator?](#)

[How much ventilation is needed for the unit after H₂O₂ decontamination and before restarting cultures?](#)

[I am reluctant to put my hand into the unit to put inner parts back after H₂O₂ decontamination. May this compromise the decontaminated condition?](#)

[Is it possible to decontaminate the connecting cable and connections of the H₂O₂ generator that connects to the chamber using H₂O₂ vapor?](#)

[Why does H₂O₂ solution remain in the H₂O₂ generator after decontamination?](#)

[How do you dispose of the remaining H₂O₂ solution?](#)

[Does H₂O₂ gas leak from the door gasket?](#)

[Where will the condensation water remain after the decontamination process?](#)

[Why does the condensation that was collected in the cold spot after decontamination and resolution contain a small amount of H₂O₂? Shouldn't the condensation be just water since all H₂O₂ gas is resolved during the process?](#)

[Does H₂O₂ affect the incubator \(including gaskets\)?](#)

[Why is the chamber temperature kept at 45°C during decontamination and resolution? And when does the chamber temperature go back to normal set temperature such as 37°C?](#)

UV Section

[What do I need to do if the UV lamp burns out during the decontamination process?](#)

[What is the use of the UV lamp "automatic extension function"?](#)

[When UV intensity degrades, can the resolving phase be executed properly?](#)

[Does the extension of the UV lamp ON time affect the performance of the incubator?](#)

CO₂ Section

[Why is a zero calibration pump no longer necessary?](#)

[What is the timing of CO₂ detection with a dual wave IR Sensor inside the chamber?](#)

[Is it necessary to clean the IR sensor?](#)

Miscellaneous

[Are the following models, MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE stackable?](#)

[When should the electric lock be opened manually?](#)



MCO-230AICUVH-PE (shown open, empty)

▣ H₂O₂ Section

What do I need to do if power failure occurs during the decontamination process?

Immediately after power recovery, the UV resolving phase (90 minutes) starts. Even after the decontamination process is completed, however, a decontamination error message will be displayed. You should start the decontamination process again. During a power failure the door remains locked.

How is H₂O₂ resolved?

H₂O₂ is resolved into water and oxygen by UV irradiation to eliminate all of the H₂O₂. Without UV irradiation, only natural degradation of H₂O₂ gas concentration occurs and the condensation water inside the chamber will have an initial concentration (6%) of H₂O₂.

What marks a resolving phase completion?

When the process is completed the concentration of H₂O₂ gas is lower than 0.1 ppm (resolving time: 90 minutes). This period is preset and based on test results, not direct monitoring of actual H₂O₂ concentration levels.

During decontamination and resolving processes, is the chamber fan operating?

Yes. The fan is required to circulate the H₂O₂ vapor evenly inside the chamber during the decontamination process.

What triggers the switch from H₂O₂ generation to the resolving phase?

Once the volume of H₂O₂ solution in the generator falls below the predetermined level as monitored by a float sensor in the H₂O₂ generator, the resolving phase starts. The misting period is approximately 10 minutes and resolving period is approximately 90 minutes.

After decontamination, is it OK to leave the incubator unattended for a night or a couple of days? (Can I start the decontamination program just before going home?)

Yes, but end user must remember to clean out residual H₂O₂ before bringing cultures back in.

Can I use a commercial H₂O₂ solution?

No. Commercial H₂O₂ solutions typically contain impurities and their concentrations of H₂O₂ are often unstable in actual measurements. We do not guarantee the parts for MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE and decontamination results when using any H₂O₂ solution other than the dedicated reagent purchased from PHCbi.

When should I perform the decontamination process?

The decontamination process should be performed whenever contamination occurs inside the chamber, or as a sterilization regiment before beginning a new culture. We also recommend performing the decontamination process just after installation and a minimum of 1x per month even if contamination isn't observed.

Why do you use a 6% MCO-H2O2-PE solution?

This is the upper limit of concentration allowable before a change in classification is necessary where it would need to be handled as a general drug in Japan and USA. The MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE incubators are designed with this specific concentration for decontamination.

What temperature is needed for the storage of the H₂O₂ reagent?

Lower than 15°C. Do not freeze because it will affect concentration.

How should I store H₂O₂ reagent (MCO-H2O2-PE)?

Before use store MCO-H2O2-PE reagent in a refrigerator. Please make sure not to freeze. Some household refrigerators may have poor temperature uniformity and regulation so please keep this in mind.

What happens if MCO-H2O2-PE is frozen? Is it OK to put it in a freezer when the expiration date has been exceeded?

We have not conducted the actual measurement, but theoretically it is presumed that the amount of H₂O₂ solution is reduced once frozen.

How is H₂O₂ reagent transported?

Air freight. Please store in a refrigerator immediately upon receipt.

Is it possible to put something that I want to decontaminate into the MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE incubators and decontaminate it with H₂O₂?

We do not guarantee the decontamination effect with other non-incubator components. This is because H₂O₂ gas may not reach all areas, grooves, or corners of a different object, which would be required for decontamination.

How is decontamination affected if extra trays or half trays are added into the incubator?

Tests have been performed that demonstrate adding 1 extra tray will not cause an issue (total 5 trays). Beyond this we have no data to confirm decontamination. If the trays used in the conventional MCO series or the half trays are put in chamber for decontamination, make sure to place them so that all surfaces are exposed.

How much ventilation is needed for the unit after H₂O₂ decontamination and before restarting cultures?

It is unnecessary to perform ventilation, since the chamber air is ventilated while putting inner parts back to their original locations after decontamination. Water condensation that remains at the bottom of the unit or under the duct which may contain low concentrations of H₂O₂ should be wiped out.

I am reluctant to put my hand into the unit to put inner parts back after H₂O₂ decontamination. May this compromise the decontaminated condition?

The purpose of H₂O₂ decontamination is to decontaminate the environment of the inner chamber, and not to maintain a sterile environment. Special facilities such as clean rooms are required to maintain a sterile environment. Any incubator with a door that allows air in and out is unable to maintain a sterile environment. High temperature sterilization incubators are subject to the same conditions.

Is it possible to decontaminate the connecting cable and connections of the H₂O₂ generator that connects to the chamber using H₂O₂ vapor?

No, it is not possible to decontaminate this by H₂O₂. However, it should not be a problem because the connection is covered by a cap during the normal cultivating operation. Therefore it is important to ensure the cap is secure.

Why does H₂O₂ solution remain in the H₂O₂ generator after decontamination?

The ultrasonic generator can be damaged if energized under dry conditions, so a bit of solution remains at the end of the cycle to prevent this from occurring.

How do you dispose of the remaining H₂O₂ solution?

Please dispose in accordance with the regulation or law of each country or state. For example, the regulation in Japan calls for disposal of H₂O₂ after diluting it with a large volume of water. After usage of the H₂O₂ equipment, the insides of the H₂O₂ generator needs to be rinsed with distilled water.

Does H₂O₂ gas leak from the door gasket?

No. If the vapor is created by heating, the air inside the chamber may leak to the outside because of volumetric expansion. However, there is no volumetric expansion for this product since the ultrasonic generator creates a cool mist.

Where will the condensation water remain after the decontamination process?

Condensation water will remain at the bottom of the duct and humidifying pan area (cold-spot).

Why does the condensation that was collected in the cold spot after decontamination and resolution contain a small amount of H₂O₂? Shouldn't the condensation be just water since all H₂O₂ gas is resolved during the process?

The condensation itself does not contain H₂O₂. However, a small amount of 6% H₂O₂ solution may splatter from the H₂O₂ generator when it is generating H₂O₂ mist, which remains as liquid and unresolved by UV. This H₂O₂ residue is mixed with the condensation, resulting in condensation containing H₂O₂ of approximately 1%.

Does H₂O₂ affect the incubator (including gaskets)?

No. The interior components are designed to handle H₂O₂ exposure.

Why is the chamber temperature kept at 45°C during decontamination and resolution? And when does the chamber temperature go back to normal set temperature such as 37°C?

Keeping the unit at 45°C improves the effectiveness of vaporizing H₂O₂ and avoids condensation on most of the interior surfaces of the chamber. Once the decontamination cycle completes, the unit will go back to the normal operating temperature.

UV Section

What do I need to do if the UV lamp burns out during the decontamination process?

Although very unlikely to occur, when the UV lamp burns out during decontamination, the electric lock remains engaged and the door can't be opened until the time frame elapses. Over time the H₂O₂ vapor will naturally dissipate to below 0.1 ppm. In this case, you must exercise caution as water condensation remains inside chamber and can include approximately 6% H₂O₂ solution. Also, DO NOT manually open the door using a key until the electrical lock is released automatically.

What is the use of the UV lamp "automatic extension function"?


To compensate for the drop in UV wavelength intensity output due to aging of the UV lamp, the incubator automatically extends the UV lamp ON time depending on the accumulated illuminated age of the bulb. The extended time is displayed as a percentage under UV timer Ext, in the UV Setting Screen. The instruction manual contains this description.

When UV intensity degrades, can the resolving phase be executed properly?

When a UV lamp life indicator blinks that means the UV lamp has been illuminated for longer than 1000 hours or 5000 hours, depending on the model of the incubator and the decontamination process cannot be started. When the lamp has less than 1000/5000 hours and the UV lamp indicator is not blinking, the lamp can resolve the H₂O₂ gas.

Does the extension of the UV lamp ON time affect the performance of the incubator?

It does not affect the performance and does not cause an increase in chamber temperature.

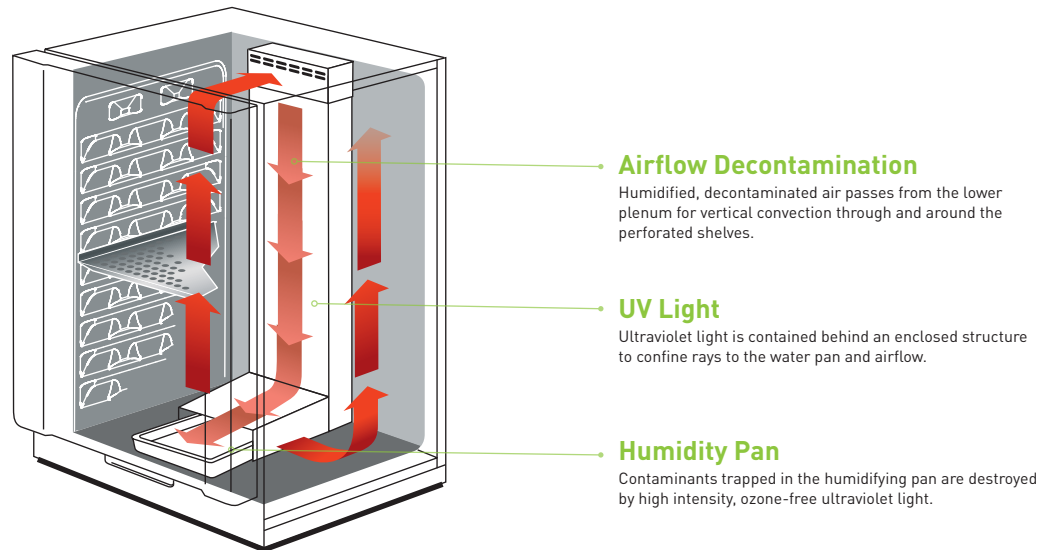

CO₂ Section

Why is a zero calibration pump no longer necessary?

The new dual wavelength IR sensor can detect CO₂ concentration inside the chamber and objective wavelength (atmosphere air as 0% CO₂) simultaneously using infrared light. This also contributes to increased reliability and lower vibration as there is no longer a need for an air pump.

What is the timing of CO₂ detection with a dual wave IR Sensor inside the chamber?

Real-time immediate response not affected by chamber humidity.

Is it necessary to clean the IR sensor?

No.

Miscellaneous

Are the following models, MCO-170AICUVH-PE, MCO-230AICUVH-PE and MCO-170MUVH-PE stackable?

Yes. See chart below. Contact PHCbi for information on stacking adapters.

Incubator Stacking Matrix for Model MCO-170AICUVH-PE / MCO-170MUVH-PE				
	Model	Model	Model	Model
Top Incubator	MCO-170AICUVH-PE	MCO-170MUVH-PE	MCO-170AICUVH-PE	MCO-170MUVH-PE
Bottom Incubator	MCO-170AICUVH-PE	MCO-170AICUVH-PE	MCO-170MUVH-PE	MCO-170MUVH-PE

Incubator Stacking Matrix for Model MCO-230AICUVH-PE			
	Model	Model	Model
Top Incubator	MCO-230AICUVH-PE	MCO-170AICUVH-PE	MCO-170MUVH-PE
Bottom Incubator	MCO-230AICUVH-PE	MCO-230AICUVH-PE	MCO-230AICUVH-PE

Even most earlier PHCbi and Sanyo incubators can be stacked on these models. Contact PHCbi for complete information.

When should the electric lock be opened manually?

The lock should be opened manually whenever the incubator is moved. The door is locked when the incubator is switched off or during a power failure. Do not open manually during decontamination operation even if a power failure occurs.

PHCbi

PHC Europe B.V.
Nijverheidsweg 120 | 4879 AZ Etten-Leur | Netherlands
T: +31 (0) 76 543 3839 | F: +31 (0) 76 541 3732
www.phchd.com/eu/biomedical